Alsident® System 100

Aluminium and Chemical Resistant
Aluminium und Chemisch Beständig
Aluminium og Kemisk Resistent

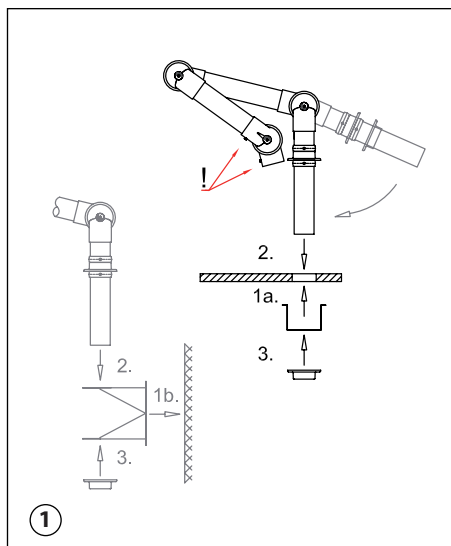
Fig. 1

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UK – System 100 TM

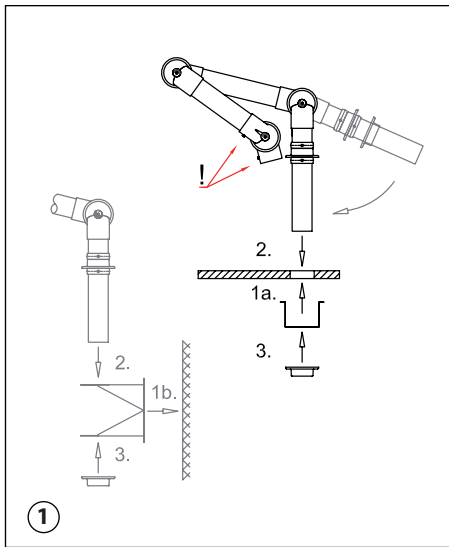
Fig. 1

1. Make a hole with a diameter of 120 mm where the extraction arm is to be placed. The U-profile is fastened to the table (1a). Or fasten the wall bracket on the wall in the desired height above the floor (1b).
2. Considering the internal spring it is important to turn the socket pipe in the direction of the arrow. Make sure the fingerscrews (!) are faces forward.
3. Place the socket pipe through the holes in the U-profile/wall bracket and fasten the flange with screws (2).
4. Place the other flange on the socket pipe under the table/bracket and fasten it with screws (3).

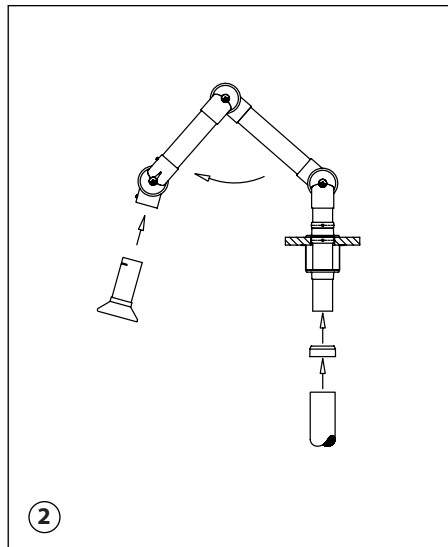
Fig. 2

1. Connect the extraction arm to the main ventilation duct with the Alsident® reducer. In order to turn the extraction arm, it is important not to fix the reducer to the socket pipe
2. Unfold the extraction arm in the direction of the arrow.
3. Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew. When screwed in the finger-screw will hold the accessory in place in the joint.
4. Finally make sure the damper is open.
5. The extraction arm is now ready for use.

Alsident® System reserve the right to make changes in this brochure to technical data, construction, materials, equipment and appearance. The illustrations do not necessarily represent the true appearance of the products and are for the purpose of description only. Alsident® System make reservations regarding misprints.

See MAINTENANCE AND CLEANING
at www.alsident.com under "Technical Information"

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